



APPLICATION NOTE

Subject : Analysis of standard material among general compounds

▶ Analysis condition

Oven : 35.0°C(5 min) → 5.0 °C/min → 150.0 °C(1min)

Injection Port 1 : Capillary; 230°C, He flow rate=1.0 mL,

Split ratio=50.0; Constant flow


Column 1 : HP-5(30m*0.32mm*1.0um)

Detector 1 : FID, 250°C

Extra condition : Diluted by DMSO. Heat on 80 °C (for 60 min.)

Injection Vol. : 1ml

▶ Chromatogram

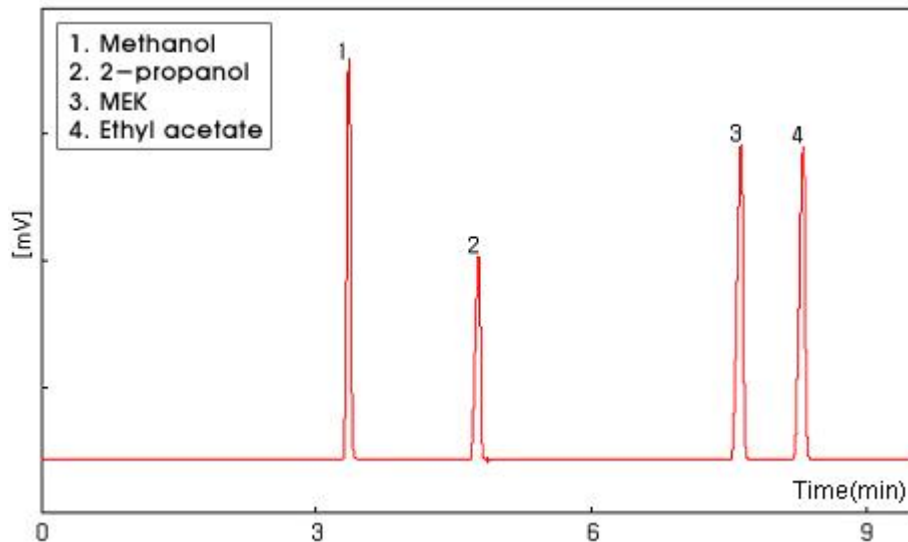

영린기기

431-836 경기도 안양시 동안구 호계2동 899-6 Tel : 031-428-8700 Fax : 031-428-8787
 Homepage : www.younglin.com E-mail : sales@younglin.com