





Analysis of Carbamate Pesticides

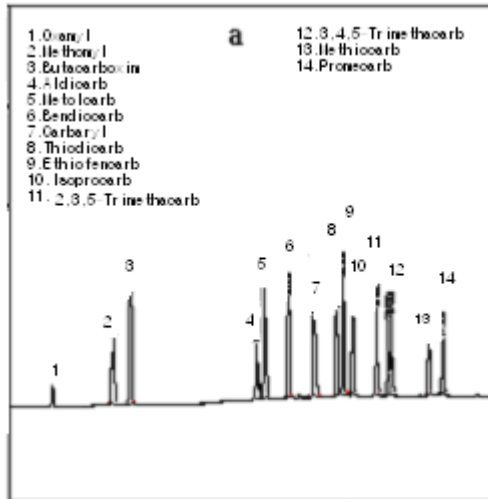

a. Carbamate Standard 5ppm

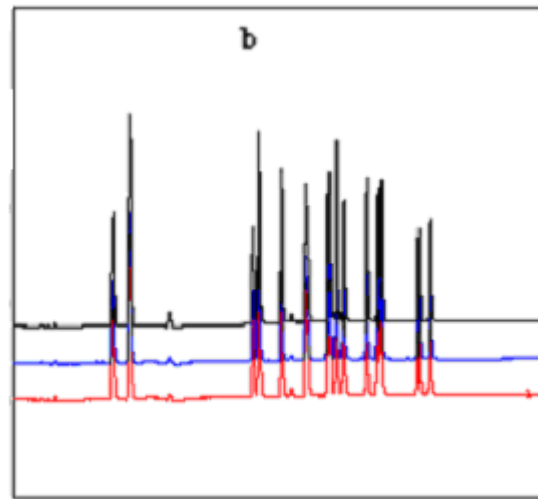

b. Overlap Chromatogram

Analysis Condition

Column : Carbamate column, 250mm X 4.6mm X 5.0um

Column temp.: 42°C

Flow rate : 1.0 mL/min

Injection vol.: 20uL

Run time: 33min

| | Water | Methanol |
|---------|-------|----------|
| Initial | 80 | 20 |
| 2 min | 70 | 30 |
| 20min | 40 | 60 |
| 25min | 30 | 70 |
| 33min | 20 | 80 |



Carbamate Analyzer