

APPLICATION NOTE

Plasma 중 카테콜아민 분석



1. 분석조건

Acme9000 Series HPLC

Mobile Phase & Column : Chromsystem kit에 포함된 제품사용


Flow Rate : 1.3ml/min

Column Oven : 35°C

Detector : ECD (+500mV, 2nA, WE: Glassy Carbon)

Injection Volumn : 20ul sample loop

2. 크로마토그램

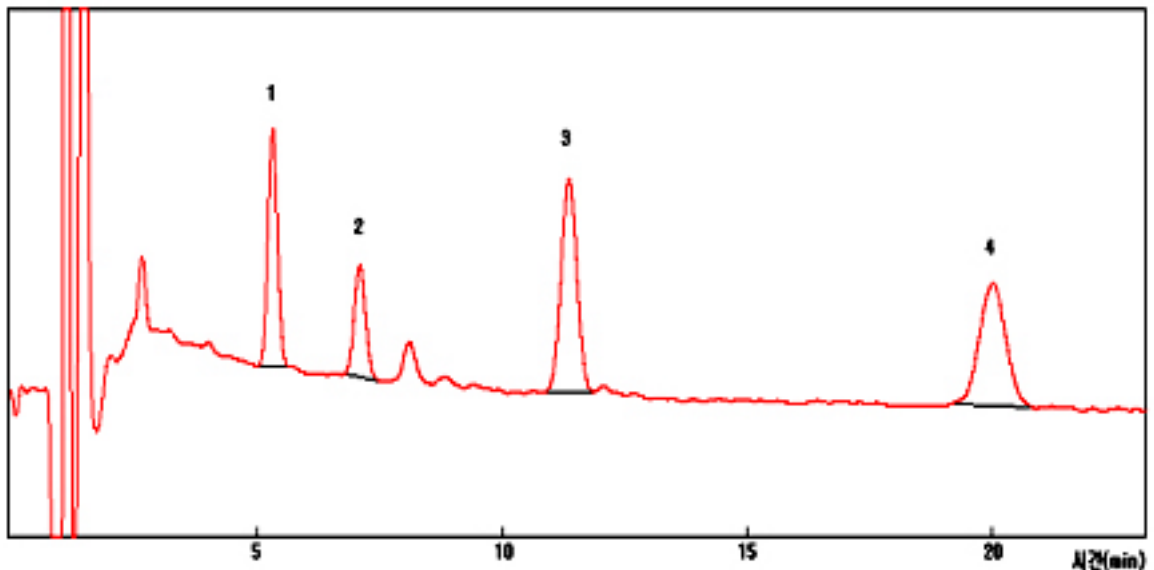

1) Norepinephrine (NE)

2) Epinephrine (Epi)

3) DHBA(ISTD)

4) Dopamine (DA)



영린기기

431-836 경기도 안양시 동안구 호계2동 899-6 Tel : 031-428-8700 Fax : 031-428-8787
 Homepage : www.younglin.com E-mail : sales@younglin.com